



pluriBead Short Protocol

universal beads

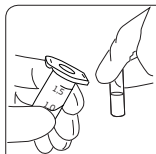


Preparation

Mix pluriBeads, your antibody and PBS solution in a 1,5ml reaction tube.

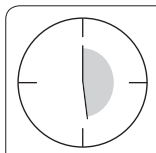


Incubation

Incubate the sample 3-4 hrs at room temperature with a thermoshaker (~750rpm). Alternatively, use mixing devices that allow for overhead tilting.

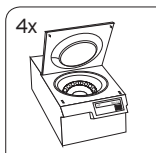

Purification

Wash with PBS, centrifuge the suspension and remove supernatant. Repeat this step 3 times.

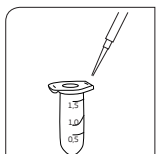

Further Usage

pluriBead universal particles are now labeled with your antibody.

Add PBS solution

See pluriBead protocol for whole blood, tissue or buffy coat for separating the required targets with your custom beads.